



## Agilent WineAnalyzer Software Installation Instructions

### NOTE

The WineAnalyzer software application is not a stand-alone program — it must be run in conjunction with at least one of the following Cary WinUV packages:

Application	Version number
Bio Pack	2.00(26) or later
Conc Pack	2.00(26) or later
Scan Pack	2.00(26) or later
Solascreen	2.00(27) or later
Tablet Pack	2.00(27) or later

It is strongly recommended that the WineAnalyzer software application be installed after the installation of one of these Cary WinUV packages.

---

### To install the WineAnalyzer software application:

- 1 If you are using the Microsoft® Windows NT® operating system, you must log on as an administrator or as a user with administrator privileges to successfully install the WineAnalyzer application software.
- 2 If you have not already done so, install the Cary WinUV software. For information on how to do this, refer to the Cary WinUV software manual (publication number 8510162500).
- 3 Insert the WineAnalyzer CD-ROM into the CD-ROM drive.
- 4 The WineAnalyzer application will automatically be installed in the :/Program Files/Agilent/ADL directory and a WineAnalyzer icon will appear on your desktop.

To run the WineAnalyzer application, double-click the WineAnalyzer icon on the desktop.



This information is subject to change without notice.



8510201400

Part Number: 8510201400

Edition 3/11

Issue 2

Printed in Australia

© Agilent Technologies, Inc. 2002, 2011

Agilent Technologies

679 Springvale Road

Mulgrave, VIC 3170