

Oligosaccharide Analysis on Agilent PLgel with Gel Permeation Chromatography

Application Note

Materials Testing and Research, Polymers

Author

Graham Cleaver
Agilent Technologies, Inc.

Introduction

Oligosaccharides are naturally occurring saccharides, typically with three to 10 component sugars. They are usually linked to lipids or amino acid side chains in proteins, and have many functions, such as cell-to-cell recognition in animals.

These compounds are soluble in water and polar organic solvents such as N-methylpyrrolidone (NMP), which may be required for more polar oligosaccharides such as starches, as this application note demonstrates. Gel permeation chromatography of oligosaccharides with NMP is straightforward using Agilent PLgel 5 μm columns.

Analysis of an Oligosaccharide

The viscosity of NMP is extremely high and operation at elevated temperature is recommended to reduce pressure and enhance resolution.



Conditions

Column	Agilent PLgel 5 μm 500Å, 7.5 \times 300 mm (p/n PL1110-6525)
Eluent	NMP
Flow rate	1.0 mL/min
Temp	80 °C
Detector	RI

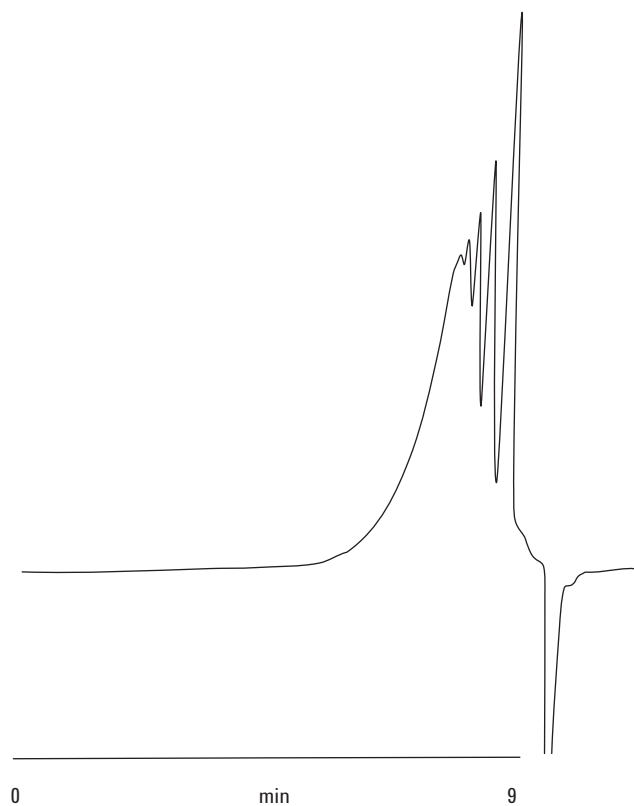


Figure 1. Separation of oligosaccharide on an Agilent PLgel 5 μm column.

Conclusions

Agilent PLgel columns in a polar solvent such as N-methylpyrrolidone can be used to analyze oligosaccharides by gel permeation chromatography.

For More Information

These data represent typical results. For more information on our products and services, visit our Web site at www.agilent.com/chem.

www.agilent.com/chem

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc., 2015
Printed in the USA
April 30, 2015
5990-8489EN