

AGILENT 7900 ICP-MS

RAISE YOUR EXPECTATIONS WITH THE NEXT GENERATION ICP-MS


Innovative technology and a new MassHunter software platform combine to make the Agilent 7900 the world's most powerful, and most automated quadrupole ICP-MS.

www.agilent.com/chem/7900icpms




WORLD'S MOST POWERFUL ICP-MS SOFTWARE

With automated system optimization, intelligent method setup, remote monitoring capability and a simplified user interface, Agilent's ICP-MS MassHunter software blends powerful features with unprecedented ease of use.

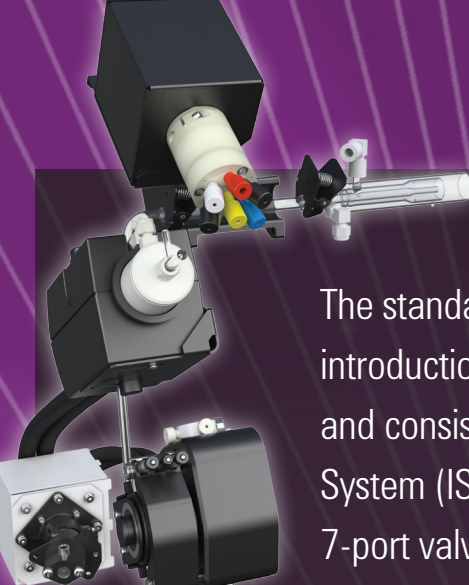

SAMPLE INTRODUCTION

The standard low-flow, Peltier-cooled sample introduction system increases operational stability and consistency. The Integrated Sample Introduction System (ISIS 3) adds a piston pump and close-coupled 7-port valve for high-speed discrete sampling.

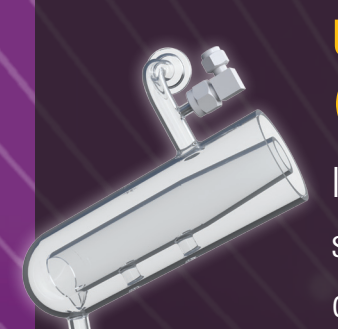

ULTRA HIGH MATRIX INTRODUCTION (UHMI)

Increases matrix tolerance up to 25% total dissolved solids (TDS). UHMI also improves plasma robustness, dramatically reducing matrix suppression.

PLASMA AND SHIELD TORCH SYSTEM (STS)

Provides precise ion energy control, ensuring high sensitivity and effective interference removal in helium mode. The torch auto-aligns with the interface following maintenance.

INTERFACE AND CONES

Standard Ni, or optional Pt-tipped cones increase ion transmission and matrix tolerance. Screw-threaded for easy removal during maintenance.




OFF-AXIS ION LENS

Improves ion transmission across the mass range, *without* the need for mass-specific voltage optimization.

4TH-GENERATION OCTOPOLE REACTION SYSTEM (ORS⁴)

Temperature-controlled collision/reaction cell with a new gas controller for fast cell gas switching in less than 3 seconds.

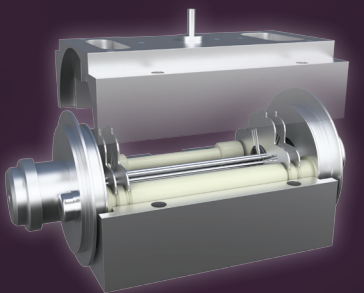

OCTOPOLE ION GUIDE


The octopole provides superior interference removal by KED in helium collision mode, and has been field-proven in thousands of Agilent ICP-MS installations.

HYPERBOLIC QUADRUPOLE

The *only* hyperbolic quadrupole used in ICP-MS. It delivers superior peak separation and abundance sensitivity, without the need for custom quadrupole settings to separate adjacent peaks.

ORTHOGONAL DETECTOR SYSTEM (ODS)

The ODS delivers higher sensitivity, lower background, and a wider measurement range — up to 11 orders of magnitude from 0.1 cps to 10 Gcps — virtually eliminating over-range results.

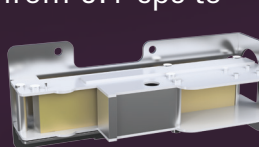

COMPACT BENCHTOP DESIGN

The world's smallest ICP-MS system saves valuable bench space while ensuring easy access for servicing and maintenance.

27 MHz PLASMA RF GENERATOR

The fast, frequency-matching RF generator increases the tolerance to changing matrices. Even volatile organic solvents can be introduced without affecting plasma stability.

VACUUM SYSTEM

A single, high-performance split-flow turbo pump and external rotary pump optimize vacuum in the interface region, increasing sensitivity while improving matrix tolerance.

AGILENT PARTS AND SUPPLIES

Manufactured to stringent specifications to ensure top quality, and rigorously tested to maximize instrument performance.



This information is subject to change without notice.

© Agilent Technologies, Inc. 2014
Printed in the USA April 1, 2014
5991-4197EN

**LEADING THE WAY IN
ATOMIC SPECTROSCOPY
INNOVATION**

www.agilent.com/chem/AtomicSpec



AGILENT ICP-QQQ




AGILENT ICP-MS



AGILENT ICP-OES



AGILENT MP-AES

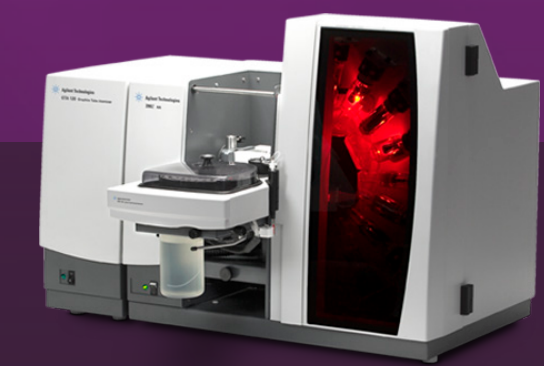

AGILENT AA

