

## Environmental analysis

# Analysis of multiple matrices with a single calibration curve for polycyclic aromatic hydrocarbons (PAHs) with the ISQ 7610 GC-MS system following EPA Method 8270E

## Authors

Chiara Calaprice<sup>1</sup>, Brian Pike<sup>2</sup>,  
Giulia Riccardino<sup>3</sup>, Adam Ladak<sup>4</sup>,  
and Paul Silcock<sup>4</sup>

<sup>1</sup>Thermo Fisher Scientific, Bremen, DE

<sup>2</sup>PACE Analytical IDEA Laboratory,  
Minneapolis, MN, USA

<sup>3</sup>Thermo Fisher Scientific, Milan, IT

<sup>4</sup>Thermo Fisher Scientific, Hemel  
Hempstead, UK

## Keywords

Polycyclic aromatic hydrocarbons (PAHs), EPA Method 8270E, ISQ 7610 mass spectrometer, extended dynamic range, gas chromatography-mass spectrometry (GC-MS), ExtractaBrite, Chromeleon, Environmental Analysis Pack

## Goal

The aim of this application note is to demonstrate the wide dynamic range and the robustness of the Thermo Scientific™ ISQ™ 7610 single quadrupole mass spectrometer, using the new Thermo Scientific™ XLXR™ detector, coupled to a Thermo Scientific™ TRACE™ 1610 gas chromatograph, for the analysis of 19 polycyclic aromatic hydrocarbons (PAHs) in soil and water, according to the United States Environmental Protection Agency (EPA) Method 8270E.

## Introduction

Polycyclic aromatic hydrocarbons (PAHs) are organic compounds consisting of carbon and hydrogen atoms. Chemically the PAHs comprise two or more aromatic rings bonded in linear, cluster, or angular arrangements, resulting in a wide diversity of physical, chemical, and toxicological properties. PAHs are ubiquitous and can contaminate soil, air, sediments, and water and are resistant to environmental degradation. These compounds are found in fossil fuel sources and manmade chemicals and are derived from the incomplete combustion of organic matter used for human activities (such as vehicle emissions, rubber, plastics, and cigarettes). PAHs have toxic effects because of their chemical structure and act as a carcinogen or endocrine disrupter. Due to their toxicity, they are monitored in the environment with strict regulations.<sup>1</sup>

One of the most common regulations followed for the analysis of PAHs is EPA Method 8270E.<sup>2</sup> Analytical laboratories following this method face several challenges. The first challenge is that isobaric compounds must have sufficient chromatographic resolution, in particular benzo[*b*]fluoranthene and benzo[*k*]fluoranthene. High boiling compounds, such as benzo[*g,h,i*]perylene, also pose a challenge as there is a possibility for carryover and peak broadening.<sup>3</sup> Careful optimization of instrumental conditions must be done to avoid saturation and linearity loss; labs may also need to separate calibration curves for different matrices, for example soil and water, to ensure they do not exceed the linear dynamic range of the system.

Following the regulations for EPA Method 8270E comes with its own challenges. DFTPP tuning must be performed to ensure that the ion abundances are acceptable for the analysis. 8270E requires a tune during the initial full calibration, then the continuing calibration to be run every 12 hours after that for analysis. All previous versions of 8270 before E required a full DFTPP tune every 12 hours. If DFTPP tune fails, the entire batch of samples must be rerun to be compliant with the method. The final challenge for analytical testing laboratories performing this analysis is to maintain the sample throughput. It is essential that the instrument performs consistently throughout the analysis, and extended runs without maintenance are desirable. If there is any unproductive time on the instrument caused by venting to clean the system or changing the column, the sampler turnaround time and asset utilization is affected and results to clients are delayed.

In this application note, the ISQ 7610 single quadrupole GC-MS system was utilized for the simultaneous analysis of PAHs in water and soil samples. The XLXR detector comes as standard on the system and offers extended linear dynamic range and lifetime. For this analysis, a single calibration curve over five orders of magnitude was utilized to analyze water and soil samples. This extended dynamic range eliminates the need to run separate curves for different matrices and aids to increase sample throughput. An extended run of soil and water matrices were also analyzed on the system to demonstrate the robustness for the analysis of PAHs. The NeverVent™ technology on the ISQ 7610 GC-MS also allows for instrument downtime to be significantly reduced due to the ability to exchange the column and clean the ionization source without needing to vent the system. By eliminating unproductive time on the instrument, more injections can be performed on the system.

## Experimental

### Reagents and standards

Native compounds calibration mix containing 18 PAHs listed in the EPA Method 8270 (each component at 2,000 µg/mL,

P/N 31995), labeled internal standard mix (2,000 µg/mL, P/N 31206) and GC-MS Tuning mix (1,000 ng/mL, P/N 31615) were purchased from Restek; neat dibenzofuran (10 mg, P/N DRE-C20710000) was obtained from LGC and diluted in dichloromethane (DCM) to a concentration of 10 mg/mL. Surrogate standard mix (4,000 µg/mL, P/N M-8270-SS) was purchased from AccuStandards.

### Preparation of solvent calibration curve, instrument detection limit (IDL), and method detection limit (MDL) samples

Thirteen calibration solutions in DCM, containing 19 native PAHs, labeled internal standards and the surrogate standard were prepared, ranging from 2.5 to 20,000 ng/mL (ppb) (full details in Appendix); the ISTD was at 1,000 ng/mL and the surrogate was at 800 ng/mL. Average response factor calibration was used, and 15% RSD criterion was applied to assess linearity in this wide calibration range. Instrument detection limit (IDL) was calculated by injecting a 2.5 ng/mL calibration solution in DCM. Method detection limit (MDL) was calculated using extracts of water and soil spiked at 2.5 and 5 ng/mL, respectively, after the extraction.

### Preparation of samples and QCs

Water and soil extracts (n=76) were provided by PACE Analytical®, USA. Samples were spiked, extracted following the EPA Methods 3510<sup>4</sup> and 3511<sup>5</sup> for water, and EPA Method 3546<sup>6</sup> for soil. Water and soil samples with low levels of PAHs were spiked to have QCs at low (0.01 ppm), middle (1 ppm), and high (10 ppm) level to check for method accuracy and robustness. All the samples were injected randomly and used to assess instrument robustness over n=150 matrix injections without inlet, column, mass spectrometer maintenance, or re-tuning.

### GC-MS/MS analysis

Liquid injections of the sample extracts were performed using a Thermo Scientific™ TriPlus™ RSH SMART autosampler. Chromatographic separation was achieved using a Thermo Scientific™ TraceGOLD™ TG-PAH 30 m × 0.25 mm i.d. × 0.10 µm column. This column allowed compliance to EPA Method 8270 in terms of resolution, as well as excellent peak shape for all the compounds, including the ones with high boiling point, due to the film thickness and the high working temperature (up to 360 °C).

For the analysis, the ISQ 7610 single quadrupole GC-MS, coupled with a TRACE 1610 GC gas chromatograph and equipped with the ExtractaBrite™ ion source, was used. The method conditions are shown in Table 1. The system was operated in Selected Ion Monitoring (SIM) mode to monitor the PAHs, and in full scan (from 50 to 500 *m/z*, dwell time 0.20 s) for the tuning solution.

**Table 1. GC-MS acquisition method parameters for the determination of 19 PAHs in water and soil samples**

Injection parameters	
Inlet module and mode	SSL, split
Liner	<a href="#">P/N 453A1925-UI</a>
Liner type and size	Thermo Scientific™ LinerGOLD™, 4 mm i.d. × 78.5 mm
Injection volume (µL)	1
Inlet temperature (°C)	300
Split flow (mL/min)	15
Carrier gas, carrier flow (mL/min), carrier mode	He, 1.5, constant flow
Split ratio	10:1
Purge flow (mL/min)	5
Pre-injection needle wash	5 times, with DCM
Post-injection needle wash	10 times with DCM, 10 times with MeOH
Chromatographic column	
Thermo Scientific™ TraceGOLD™ TG-PAH	<a href="#">P/N 26055-0470</a>
Column dimensions	30 m × 0.25 mm i.d. × 0.10 µm
Oven temperature program	
Temperature 1 (°C)	40
Hold time (min)	1
Temperature 2 (°C)	285
Rate (°C/min)	35
Temperature 3 (°C)	295
Rate (°C/min)	3
Temperature 4 (°C)	350
Rate (°C/min)	30
Hold time (min)	2
Total GC run time (min):	15.2
MS parameters	
Ion source	ExtractaBrite
Transfer line temperature (°C)	350
Ion source temperature (°C)	350
Ionization type	EI
Electron energy (eV)	70
Emission current (µA)	10
Acquisition mode	SIM, 2 ions/compound
Dwell time (s)	0.02

A full list of analytes, as well as quantifier and qualifier ions that were monitored, is listed Table 2. The system was tuned with a built-in EPA Method 8270E specifically designed tune type.

**Table 2. List of target PAH compounds, with their retention time and SIM quantification and confirmatory ions**

Compound name	Rt (min)	MS quantifier ion (m/z)	MS confirmatory ion (m/z)
Naphthalene-d <sub>8</sub>	4.7	136	108
Naphthalene	4.8	128	129
2 - methyl Naphthalene	5.2	142	141
1 - methyl Naphthalene	5.3	142	141
Acenaphthylene	5.9	152	151
Acenaphthene	6.0	153	154
Acenaphthene-d <sub>10</sub>	6.0	162	164
Dibenzofuran	6.1	168	139
Fluorene	6.4	165	166
Phenanthrene-d <sub>10</sub>	7.2	188	184
Phenanthrene	7.2	178	176
Anthracene	7.2	178	176
Fluoranthene	8.1	202	200
Terphenyl-d <sub>14</sub>	8.3	244	122
Pyrene	8.4	202	200
Benz[a]anthracene	9.5	228	226
Chrysene-d <sub>12</sub>	9.7	240	236
Chrysene	9.7	228	226
Benzo[b]fluoranthene	11.3	252	250
Benzo[k]fluoranthene	11.4	252	250
Benzo[a]pyrene	12.1	252	250
Perylene-d <sub>12</sub>	12.2	264	260
Dibenzo[a,h]anthracene	13.5	278	139
Indeno[1,2,3-cd]pyrene	13.5	276	138
Benzo[g,h,i]perylene	13.9	276	138

### Data processing

Data were acquired, processed, and reported using Thermo Scientific™ Chromeleon™ 7.3 CDS software, which allows instrument control, method development, quantitative/qualitative analysis, and customizable reporting all within one platform. The GC-MS Environmental Extension Pack includes a suite of report templates, processing methods and eWorkflows™ to facilitate environmental analysis by GC-MS using EPA Methods 8270,

524, 525, and 8260<sup>6</sup>. For the analysis of PAHs, monitoring the ratios between the masses of the DFTPP is required and automatically performed by the software System Suitability Check. Chromeleon CDS allows rapid implementation of the PAHs method into any analytical laboratory and ensures the system produces results shortly after installation.


## Results and discussion

### Chromatographic separation and resolution of isomers

The optimized GC conditions and the high selectivity of the TraceGOLD TG-PAH capillary column allowed for chromatographic resolution of isobaric compounds in a total run time of 15.2 minutes, meeting the EPA Method 8270E requirements. Gaussian peak shapes were obtained for all the compounds, including the ones with high boiling points.


An example chromatogram obtained for a solvent standard at 0.1 ppm is shown in Figures 1 and 2. Peaks of naphthalene and benzo[*g,h,i*]pyrene, the first and the last eluting PAHs, show gaussian and sharp peak shape from the beginning to the end of the method; baseline peak width for benzo[*g,h,i*]pyrene was 0.034 min.

Phenanthrene and anthracene compounds are almost baseline resolved; benzo[*b*]fluoranthene and benzo[*k*]fluoranthene have a resolution of 20% (calculated as the ratio between the height of the valley and the smaller height of the apex of the two compounds); dibenzo[*a,h*]anthracene and indeno[1,2,3-*cd*]pyrene coelute and rely on mass separation.

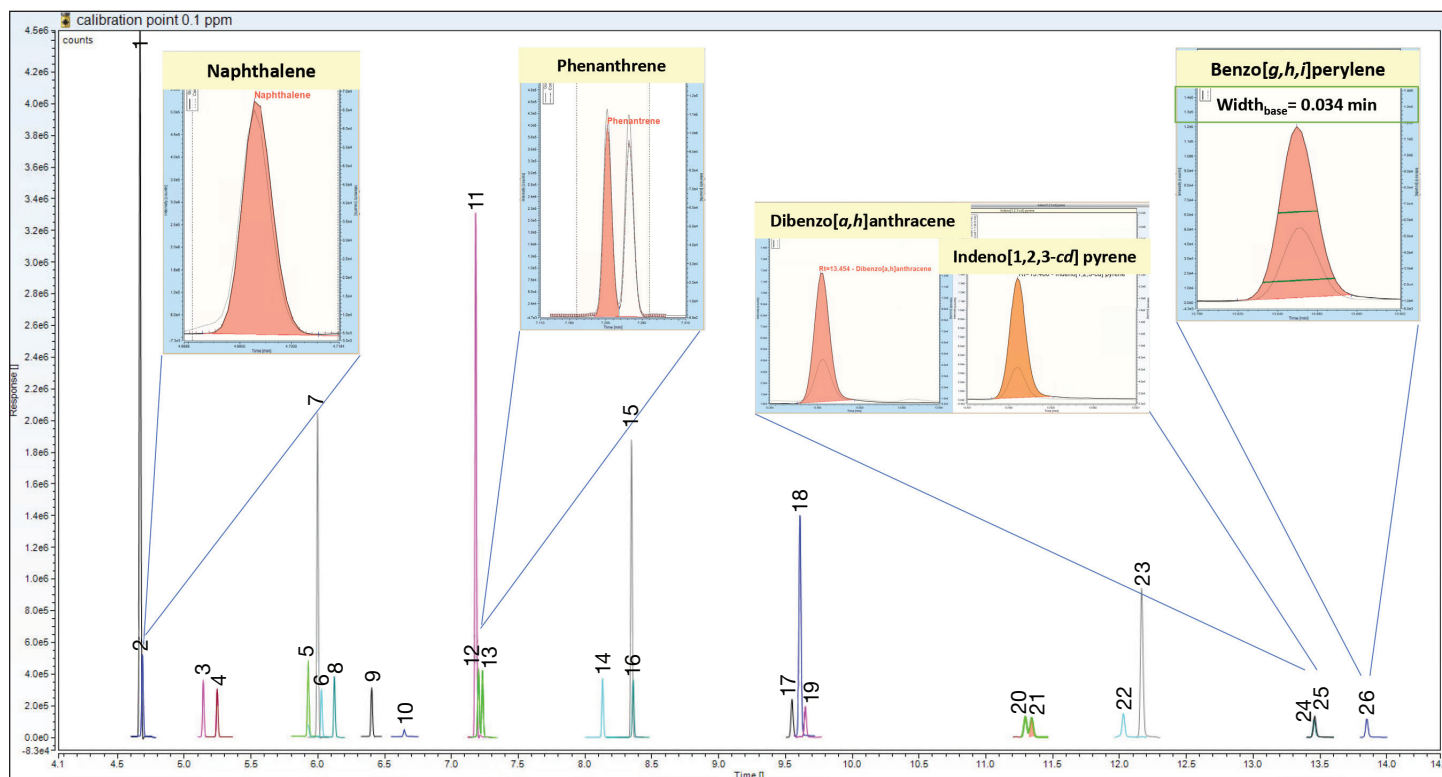

1 = Naphthalene-*d*<sub>8</sub>, 2 = Naphthalene, 3 = 2-methyl Naphthalene, 4 = 1-methyl Naphthalene, 5 = Acenaphthylene, 6 = Acenaphthene, 7 = Acenaphthene-*d*<sub>10</sub>, 8 = Dibenzofuran, 9 = Fluorene, 10 = Tribromophenol 2,4,6, 11 = Phenanthrene-*d*<sub>10</sub>, 12 = Phenanthrene, 13 = Anthracene, 14 = Fluoranthene, 15 = Terphenyl-*d*<sub>14</sub>, 16 = Pyrene, 17 = Benz[*a*]anthracene, 18 = Chrysene-*d*<sub>12</sub>, 19 = Chrysene, 20 = Benzo[*b*]fluoranthene, 21 = Benzo[*k*]fluoranthene, 22 = Benzo[*a*]pyrene, 23 = Perylene-*d*<sub>12</sub>, 24 = Dibenzo[*a,h*]anthracene, 25 = Indeno[1,2,3-*cd*]pyrene, 26 = Benzo[*g,h,i*]pyrene

Figure 1. Chromatographic separation and peak shape for 19 investigated PAHs, 5 labeled ISTD, and 2 surrogate standards in a solvent standard at 0.1 ppm acquired in SIM. Tracks of the quantification ions are reported for each peak in different colors.

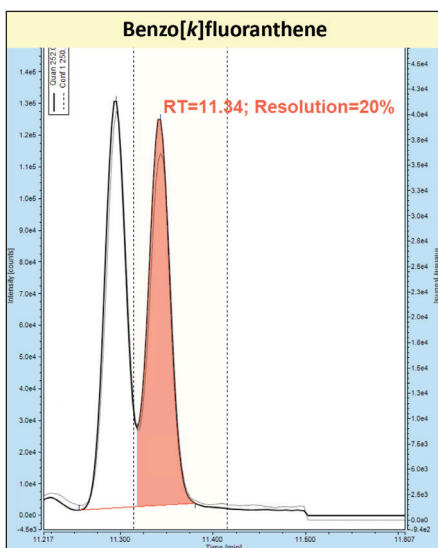


**Figure 2. Chromatographic separation of benzo[b]fluoranthene and benzo[k]fluoranthene.** Peak resolution is calculated as the ratio between the height of the peak valley and the height of the peak apex of the smaller compound.

**Table 3. Average Relative Response Factors (AvRF) calculated for each compound, and %RSD calculated over the 13 calibration points, from 2.5 to 20,000 ng/mL.** Compounds are grouped by reference ISTD.

Compound name	Category	AvRFs	RSD %
Naphthalene-d <sub>8</sub>	ISTD		
Naphthalene		1.16	6.60
2 - methyl Naphthalene		0.80	5.68
1 - methyl Naphthalene		0.69	7.74
Acenaphthylene		2.51	9.43
Acenaphthene		1.57	8.69
Acenaphthene-d <sub>10</sub>	ISTD		
Dibenzofuran		1.99	7.00
Fluorene		1.65	9.05
Phenanthrene-d <sub>10</sub>	ISTD		
Phenanthrene		1.28	6.00
Anthracene		1.25	4.88
Fluoranthene		1.21	7.10
Terphenyl-d <sub>14</sub>	Surrogate		
Pyrene*		2.17	8.78
Benz[a]anthracene		1.85	6.84
Chrysene-d <sub>12</sub>	ISTD		
Chrysene		1.48	5.64
Benzo[b]fluoranthene		1.63	4.85
Benzo[k]fluoranthene		1.55	5.12
Benzo[a]pyrene		1.70	6.06
Perylene-d <sub>12</sub>	ISTD		
Dibenzo[a,h]anthracene		1.20	5.94
Indeno[1,2,3-cd]pyrene		1.48	7.64
Benzo[g,h,i]perylene		1.25	7.01

## Calibration and linearity

A calibration curve ranging from 2.5 ng/mL to 20,000 ng/mL was prepared to assess linearity with the average relative response factor (AvRF) calibration fit type and the %RSD <15% criterion required in the EPA Method 8270E. Calculated %RSDs were <10% for all the compounds across the entire calibration range. The extended dynamic range of the XLXR detector allowed for using only one calibration curve rather than two different ones for the analysis of trace level and highly contaminated samples. Calculated AARF and %RSD results are reported in Table 3. Examples of calibration curves are shown in Figure 3.

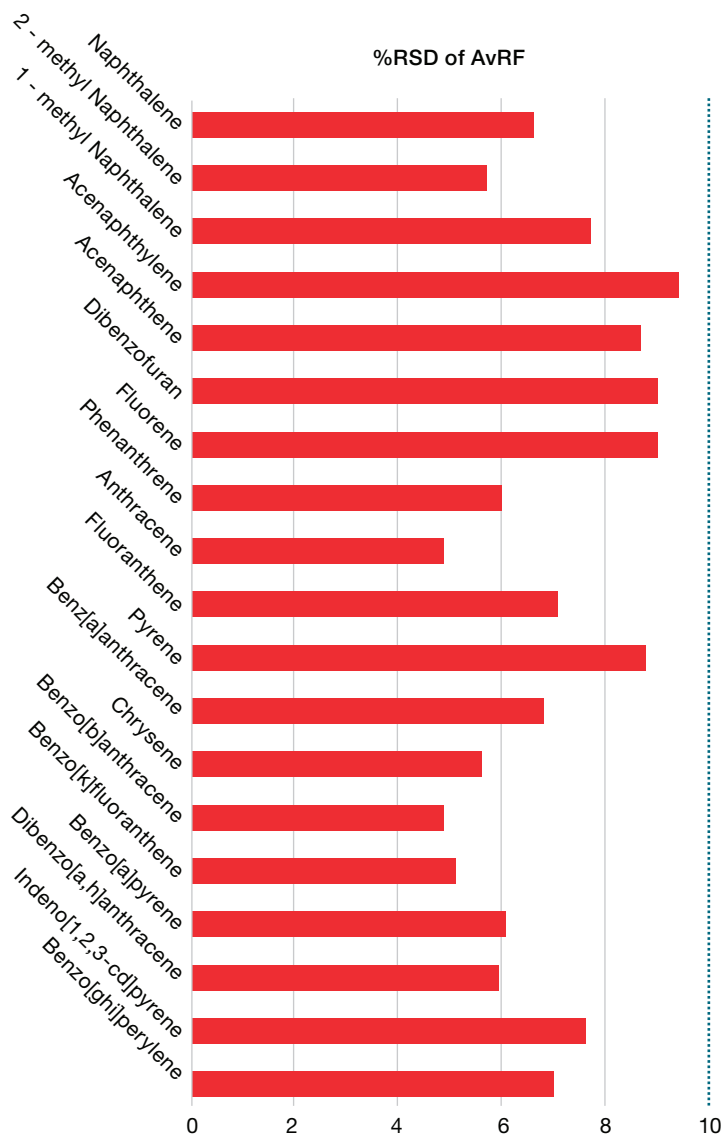

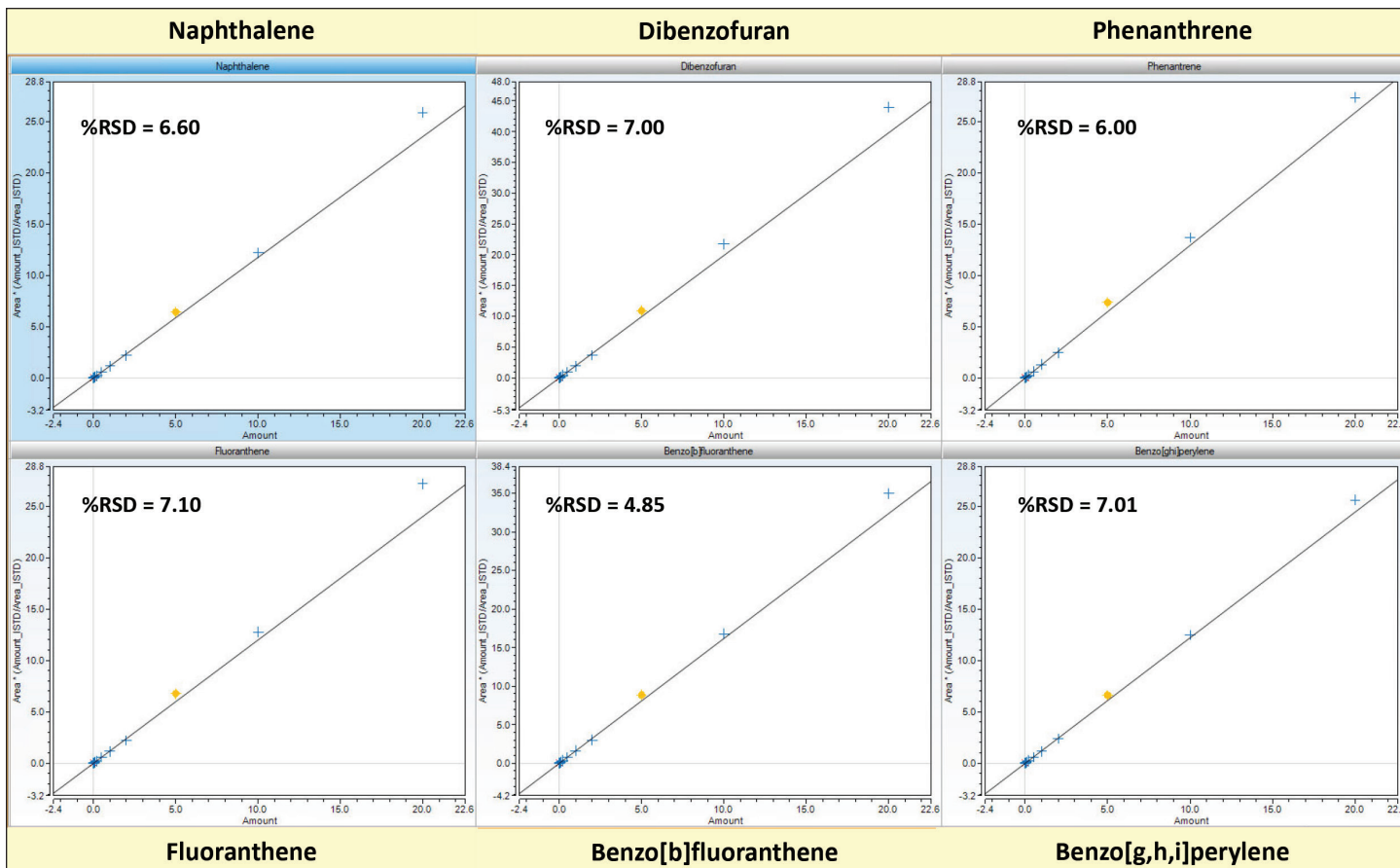


Figure 3. Examples of calibration curve for some of the 19 PAHs analyzed (from left to right: naphthalene, dibenzofuran, phenanthrene, fluoranthene, benzo[b]fluoranthene and benzo[g,h,i]perylene), annotated with the Average Response Factor %RSD. Linearity spans over five orders of magnitudes, from 2.5 to 20,000 ng/mL.

### Sensitivity: Instrument detection limits (IDLs) and method detection limit (MDLs).

To determine the instrument detection limit (IDL) for each analyte, the response and its standard deviation were taken into account.<sup>8</sup> This approach gives much more reliable LOD values as it considers not only the signal intensity, but also the consistency of the response. IDLs were calculated by eight repeated injections of the 2.5 ng/mL calibration solution in DCM, using Equation 1.<sup>8</sup>

$$IDL = t_{(n-1, 1-\alpha=0.99)} \times mOC \times \%RSD \quad (\text{Equation 1})$$

Where:

$t$  = Student  $t$ -value for one-tailed distribution at 99% confidence with 7 degrees of freedom (2.998)

mOC = mass of analyte on column

%RSD = %RSD of the response (peak area) over the 8 repeated injections. %RSD must be consistent (%RSD < 15%).

IDLs ranged from 0.3 to 0.97 pg on columns, as reported in Table 4.

The method detection limit (MDL) is defined as the minimum measured concentration of a substance that can be reported with 99% confidence that the measured concentration is distinguishable from method blank results.<sup>9</sup> For the calculation of MDL, blank extracts of water and soil were selected, pooled, and spiked at different levels (2.5 ng/mL and 5 ng/mL) after the extraction; eight repeated injections were made, and the calculations were done applying the formula reported in Equation 2.

$$MDL_s = t_{(n-1, 1-\alpha=0.99)} \times S_s \quad (\text{Equation 2})$$

Where:

$t$  = Student  $t$ -value for one-tailed distribution at 99% confidence with 7 degrees of freedom (2.998)

$S_s$  = sample standard deviation of the calculated concentration

MDLs ranged from 0.44 to 6.94 pg on column for water, and from 0.48 to 7.60 pg on column for soil.

**Table 4. Instrument detection limit in pg on column for the 19 native PAHs in DCM solvent**

Compound name	IDLs pg on column	MDLs water pg on column	MDLs soil pg on column
Naphthalene	0.97	0.44	0.48
2 - methyl Naphthalene	0.30	0.42	0.64
1 - methyl Naphthalene	0.42	0.71	0.51
Acenaphthylene	0.59	0.40	0.73
Acenaphthene	0.51	2.27	0.81
Dibenzofuran	0.42	0.61	0.65
Fluorene	0.58	1.04	1.46
Phenanthrene	0.45	1.69	1.33
Anthracene	0.64	2.92	1.40
Fluoranthene	0.39	1.84	1.73
Pyrene	0.58	1.57	1.94
Benz[ <i>a</i> ]anthracene	0.49	0.66	0.76
Chrysene	0.33	0.49	3.45
Benzo[ <i>b</i> ]fluoranthene	0.37	1.59	5.71
Benzo[ <i>k</i> ]fluoranthene	0.50	1.11	7.60
Benzo[ <i>a</i> ]pyrene	0.51	6.94	6.59
Dibenzo[ <i>a,h</i> ]anthracene	0.67	1.24	1.76
Indeno[1,2,3- <i>cd</i> ]pyrene	0.41	1.73	4.36
Benzo[ <i>g,h,i</i> ]perylene	0.33	0.86	1.48

## Repeatability

Repeatability was tested by n=10 consecutive injections of a 20 ng/mL QC in water and 20 ng/mL QC in soil. The QCs were prepared spiking blank matrices. Results are reported in Figure 4. Overall %RSD was <5.5% for water and <15% for soil matrix.

## Instrument robustness for the analysis of PAHs in environmental samples according to EPA Method 8270

System robustness and suitability for high-throughput analysis was assessed for a total of n=136 samples and 52 hours of consecutive injections, without any GC and MS maintenance (liner change, column trimming, MS tune). Details on the batch are reported in Figure 5.

## Tune check

Tuning verification was performed after each series of samples, every 6 hours, injecting 1 µL of a tuning verification solution containing DFTPP at 20 µg/mL, acquired in full scan. Tune verification criteria reported in EPA Method 8270 were automatically verified by the system suitability test (SST) in Chromeleon CDS (Figure A1 in Appendix). At the end of the batch, tune verification was still passing the specifications.

## Calibration check

Calibration check throughout the batch was done by injecting the calibration standards at different levels. The last two samples of the batch were the lowest (2.5 ng/mL) and the second to lowest (5 ng/mL) calibration points. (The accuracy in the back-calculation of these concentrations was still within ±10%.)

## QCs

QCs for water and soil matrices, at three different levels (10, 1,000, and 10,000 ng/mL) were prepared and injected every 17 samples. Calculated concentrations were accurate and consistent over the batch and were ±20% of the amount spiked (Figures 6 and 7).

### Repeatability

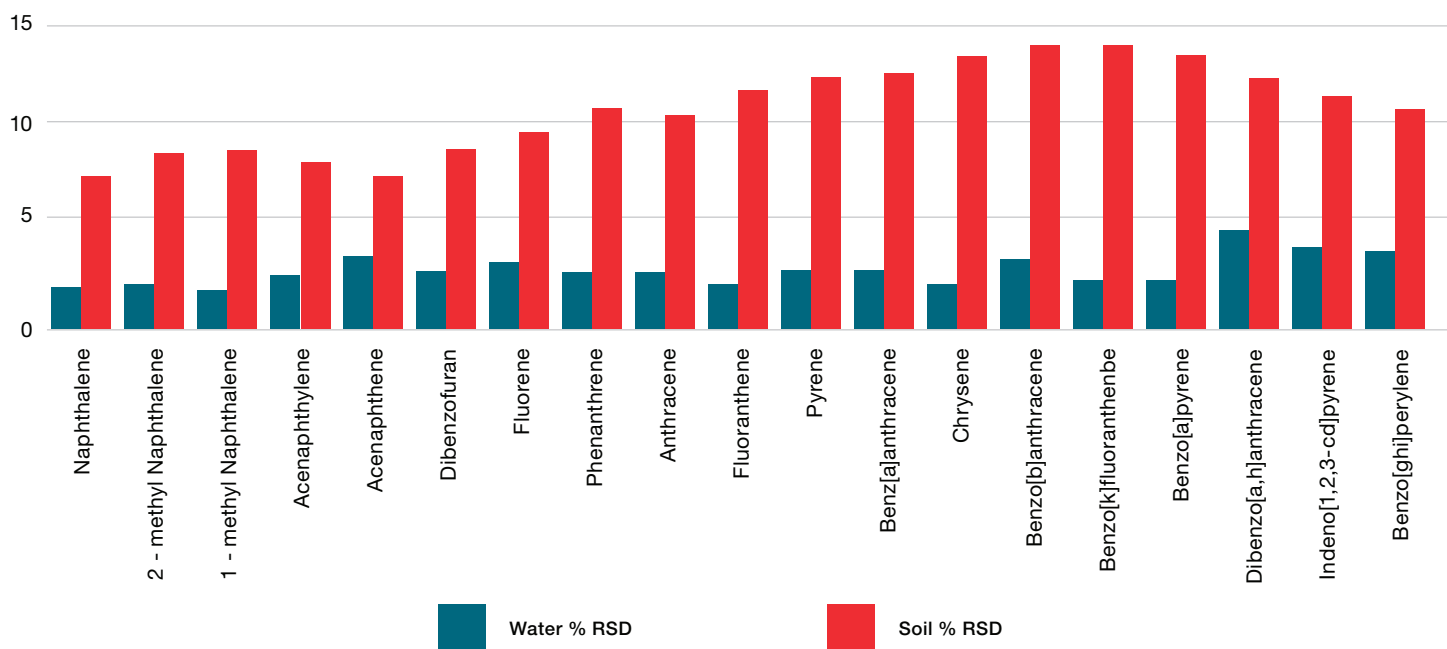


Figure 4. Peak area precision calculated in blank water and blank soil spiked at 20 ng/mL

#	TIC	Name	Type	Position	Instrument Method	Status	*Analysis_Type
1		DCM	Unknown	54	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
2		tuning mix 1	Unknown	53	PAHs SIM - 10uA - SPLIT 10to1 - FS	Finished	Field Sample
3		QC low water	Unknown	1	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
4		QC low soil	Unknown	2	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
5		QC middle water	Unknown	3	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
6		QC middle soil	Unknown	4	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
7		QC high water	Unknown	5	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
8		QC high soil	Unknown	6	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
9		Cali check 0.0025	Unknown	10	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
10		s1	Unknown	55	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
11		s2	Unknown	56	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
12		s3	Unknown	57	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
13		s4	Unknown	58	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
14		s5	Unknown	59	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
15		s6	Unknown	60	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
16		s7	Unknown	61	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
17		s8	Unknown	62	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
18		s9	Unknown	63	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
19		tuning mix 2	Unknown	53	PAHs SIM - 10uA - SPLIT 10to1 - FS	Finished	Field Sample
20		Cali check 0.005	Unknown	11	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
21		QC low water 2	Unknown	1	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
22		QC low soil 2	Unknown	2	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
23		QC middle water 2	Unknown	3	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
24		QC middle soil 2	Unknown	4	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
25		QC high water 2	Unknown	5	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
26		QC high soil 2	Unknown	6	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
27		s10	Unknown	64	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample
28		s11	Unknown	65	PAHs SIM - 10uA - SPLIT 10to1	Finished	Field Sample

Figure 5. Part of the batch used for robustness assessment. The batch contained 8 series of 17 samples each, namely 1 tuning mix, 1 calibration verification (CCV), 6 QCs (QC low, middle, and high for water and soil matrices) and 9 extracts of water and soil, randomly chosen. A total of n=136 samples were injected, for 52 hours of consecutive injections.



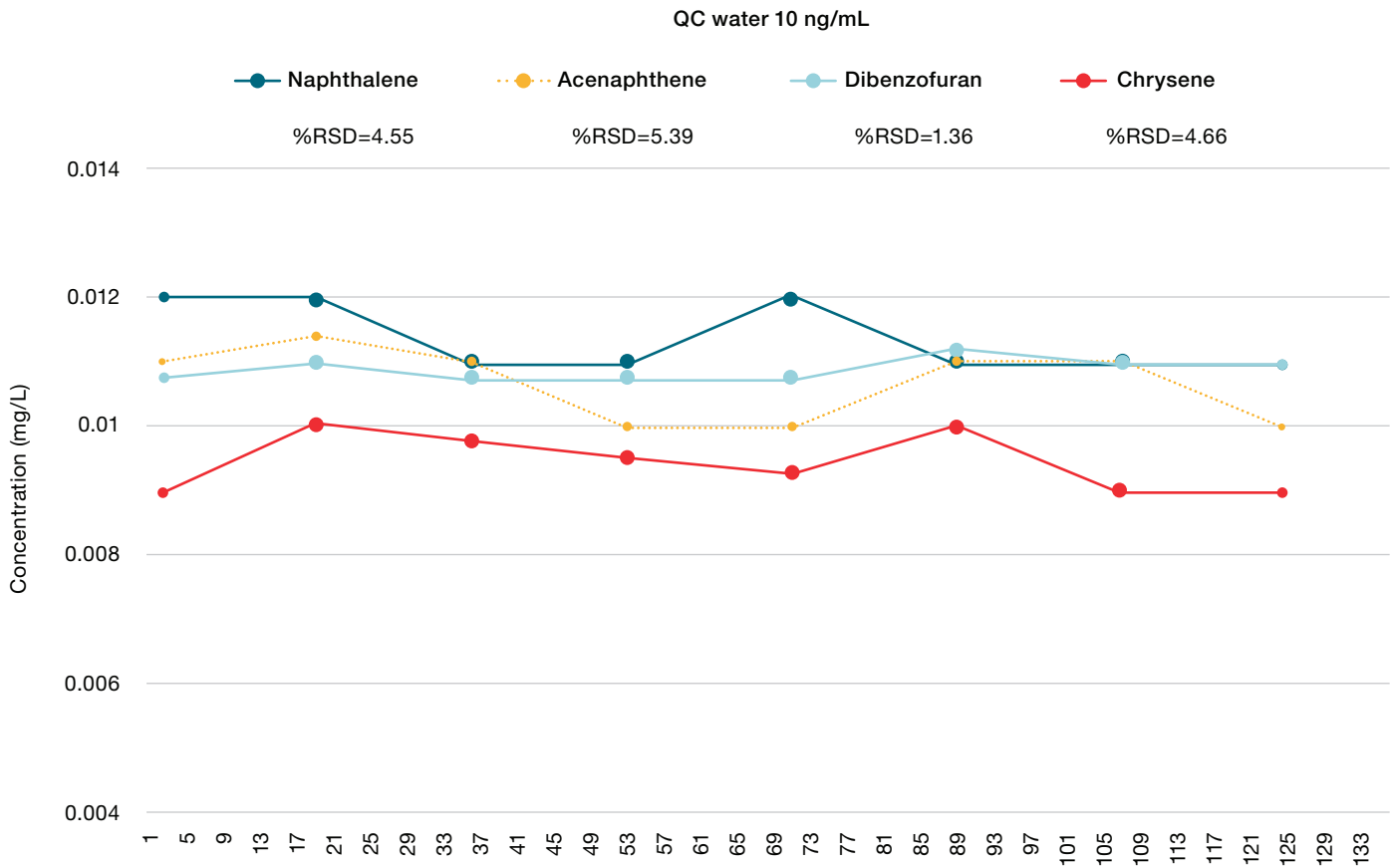


Figure 6. Calculated value of QC water blank matrix spiked at 10 ng/mL

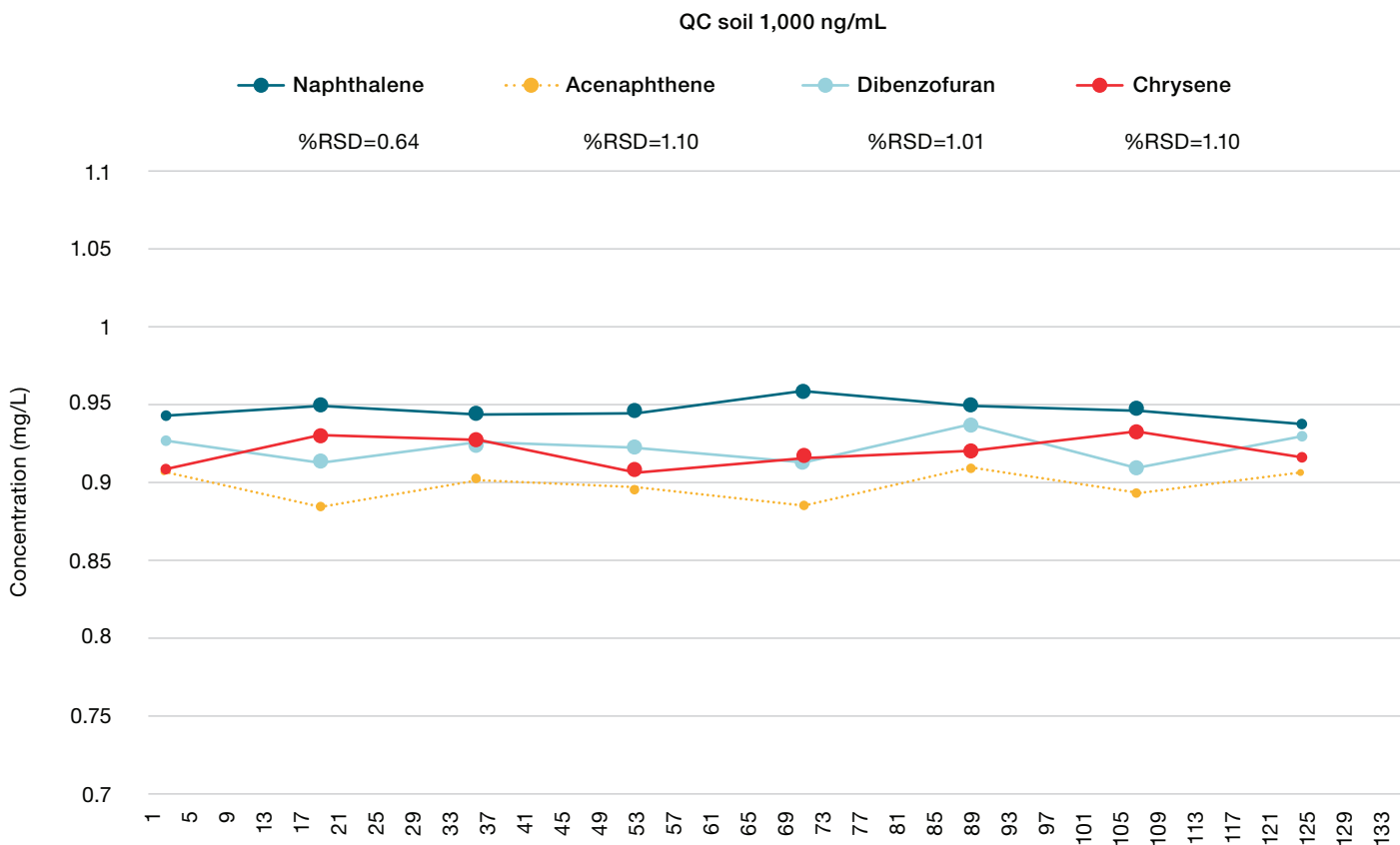


Figure 7. Calculated value of QC soil, blank matrix spiked at 1,000 ng/mL. Calculated values are consistent and accurate ( $\pm 15\%$ ) after 52 consecutive hours of injections and  $n=136$  samples.

## Conclusions

The results of the experiments described here demonstrate the ISQ 7610 GC-MS system meets and exceeds the EPA Method 8270E requirements for the analysis of PAHs in environmental samples.

- Gaussian peak shape and chromatographic resolution were achieved in compliance to EPA Method 8270E.
- The XLXR detector demonstrated exceptional dynamic range, spanning over 4 orders of magnitude, from 2.5 ppb to 20 ppm for the 19 PAHs, with AvRF %RSD<10% for each compound, thus removing the need for the preparation of two different calibration curves and speeding up the analysis workflow.
- IDLs and MDLs were in the low pg on column range, even for spiked matrices.
- The system was shown to be robust, as performances were unchanged in terms of sensitivity, peak shape, and DFTPP tuning check, after 136 injections and 52 hours of continuous work.
- The ISQ 7610 GC-MS system with NeverVent technology allows for ion source cleaning and column replacement without breaking the vacuum, thus ensuring minimal instrument downtime and high sample throughput. Unproductive instrument time is eliminated so more tests can be performed on the system.

## References

1. Lammel, G. Polycyclic Aromatic Compounds in the Atmosphere – A Review Identifying Research Needs. *Polycyclic Aromatic Compounds* **2015**, *35*, 316–329.
2. US EPA Method 8270E: Semivolatile Organic Compounds by Gas Chromatography/Mass Spectrometry, Revision 6, June 2018.
3. Thermo Scientific Application Note 10718: A fast, robust method for routine determination of polycyclic aromatic hydrocarbon (PAHs) in drinking water by single quadrupole GC-MS. Michela Balsamo *et al.* <https://assets.thermofisher.com/TFS-Assets/CMD/Application-Notes/an-10718-gc-ms-pah-drinking-water-an10718-en.pdf>
4. US EPA Method 3510: Separatory Funnel Liquid-Liquid Extraction, December 1996.
5. US EPA Method 3511 (SW-846): Organic Compounds in Water by Microextraction, Revision 1, 2014.
6. US EPA Method 3546: Microwave Extraction, February 2007.
7. Thermo Scientific Technical Note 10721: Routine, robust analysis of PAHs and PCBs in soil by single quadrupole GC-MS. Aaron Lamb *et al.* <https://assets.thermofisher.com/TFS-Assets/CMD/Technical-Notes/tn-10721-gc-ms-pahs-pcbs-soil-tn10721-en.pdf>
8. Thermo Scientific Technical Note 10499: Practical determination and validation of instrument detection limit of Thermo Fisher Scientific ISQ LT Single Quadrupole GC-MS. Richard Law, Tommaso Albertini. <https://assets.thermofisher.com/TFS-Assets/CMD/Technical-Notes/tn-10494-single-quadrupole-gc-ms-tn10494-en.pdf>
9. US EPA: Definition and Procedure for the Determination of the Method Detection Limit, Revision 2, December 2016.

# Appendix

Table A1. Concentration levels in the calibration curve

Compound name	Concentration native compounds ng/mL (ppb)	Concentration ISTDs ng/mL (ppb)	Concentration surrogate standard ng/mL (ppb)
1	2.5	1,000	800
2	5	1,000	800
3	10	1,000	800
4	20	1,000	800
5	50	1,000	800
6	100	1,000	800
7	200	1,000	800
8	500	1,000	800
9	1,000	1,000	800
10	2,000	1,000	800
11	5,000	1,000	800
12	10,000	1,000	800
13	20,000	1,000	800

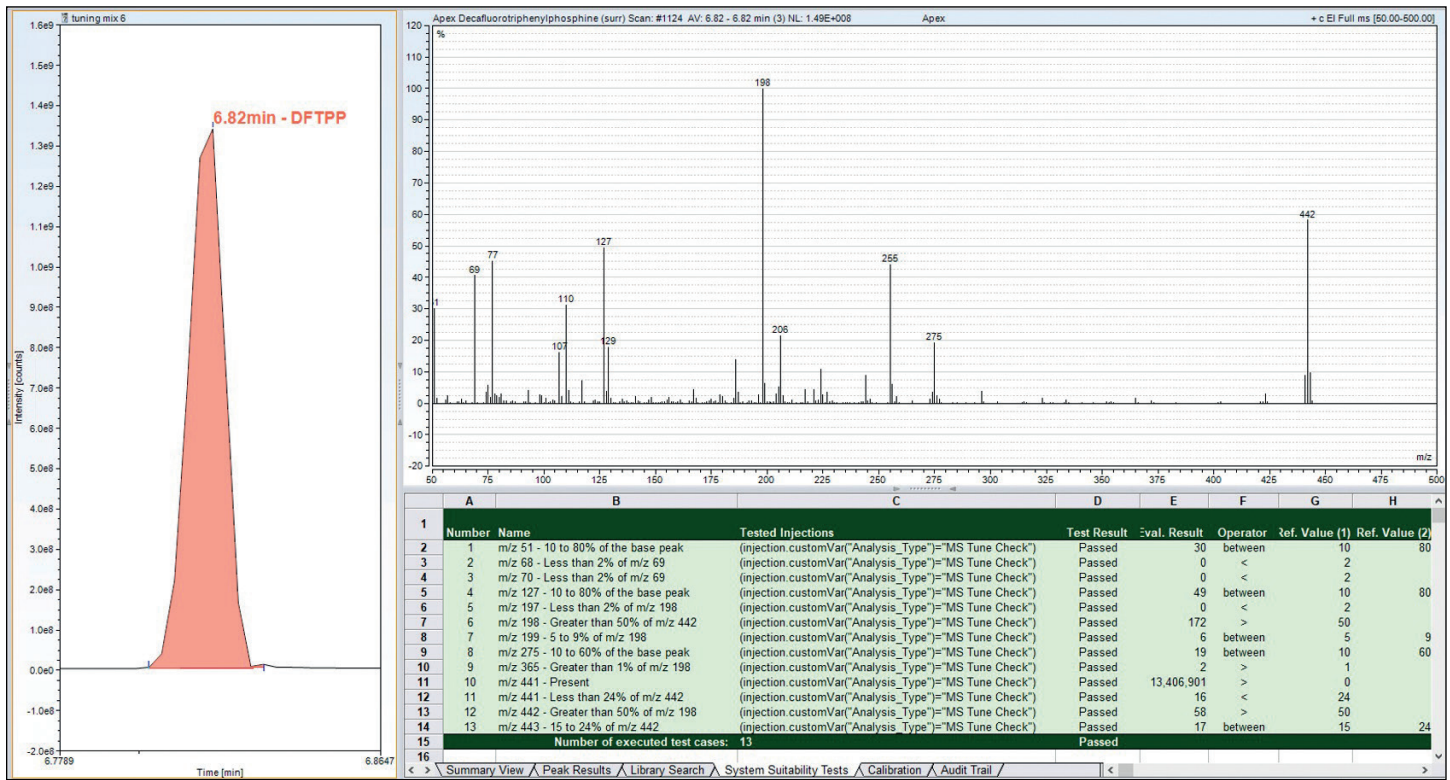


Figure A1. System suitability test results for DFTPP in the Chromeleon Data System– Environmental Pack

Table A2. Calculated concentrations of QC water at 10 ng/mL during the robustness study, accuracy, standard deviation, and %RSD

QC water 10 ng/mL	Calculated value									Target	Accuracy	2*STDEV	%RSD
Naphthalene	0.012	0.012	0.011	0.011	0.012	0.011	0.011	0.011	0.0114	0.0100	114	0.00104	4.55
2 - methyl Naphthalene	0.011	0.011	0.011	0.011	0.011	0.01	0.011	0.01	0.0108	0.0100	108	0.00093	4.31
1 - methyl Naphthalene	0.01	0.01	0.01	0.01	0.01	0.009	0.01	0.009	0.0098	0.0100	98	0.00093	4.75
Acenaphthylene	0.009	0.009	0.009	0.009	0.009	0.008	0.008	0.008	0.0086	0.0100	86	0.00104	6.00
Acenaphthene	0.011	0.0114	0.011	0.01	0.01	0.011	0.011	0.01	0.0107	0.0100	107	0.00115	5.39
Dibenzofuran	0.0108	0.011	0.0108	0.0108	0.0108	0.0112	0.011	0.011	0.0109	0.0100	109	0.00030	1.36
Fluorene	0.012	0.013	0.012	0.012	0.012	0.012	0.012	0.011	0.0120	0.0100	120	0.00107	4.45
Phenanthrene	0.038	0.037	0.038	0.037	0.037	0.038	0.036	0.036	0.0371	0.0100	371	0.00167	2.25
Anthracene	0.011	0.013	0.012	0.012	0.011	0.013	0.011	0.013	0.0120	0.0100	120	0.00185	7.72
Fluoranthene	0.02	0.019	0.02	0.02	0.019	0.019	0.019	0.019	0.0194	0.0100	194	0.00104	2.67
Pyrene	0.013	0.014	0.014	0.015	0.015	0.014	0.013	0.013	0.0139	0.0100	139	0.00167	6.01
Benzo[a]anthracene	0.009	0.009	0.009	0.009	0.009	0.009	0.01	0.01	0.0093	0.0100	93	0.00093	5.00
Chrysene	0.009	0.01	0.009	0.009	0.009	0.01	0.009	0.009	0.0093	0.0100	93	0.00093	5.00
Benzo[b]fluoranthene	0.009	0.009	0.009	0.009	0.008	0.009	0.009	0.008	0.0088	0.0100	88	0.00093	5.29
Benzo[k]fluoranthene	0.009	0.008	0.009	0.009	0.008	0.008	0.008	0.008	0.0084	0.0100	84	0.00104	6.18
Benzo[a]pyrene	0.009	0.011	0.01	0.01	0.011	0.01	0.009	0.009	0.0099	0.0100	99	0.00167	8.45
Dibenzo[a,h]anthracene	0.009	0.01	0.011	0.009	0.01	0.01	0.01	0.01	0.0099	0.0100	99	0.00128	6.49
Indeno[1,2,3-cd]pyrene	0.009	0.01	0.01	0.009	0.01	0.01	0.01	0.01	0.0098	0.0100	98	0.00093	4.75
Benzo[ghi]perylene	0.009	0.01	0.011	0.009	0.011	0.01	0.01	0.01	0.0100	0.0100	100	0.00151	7.56

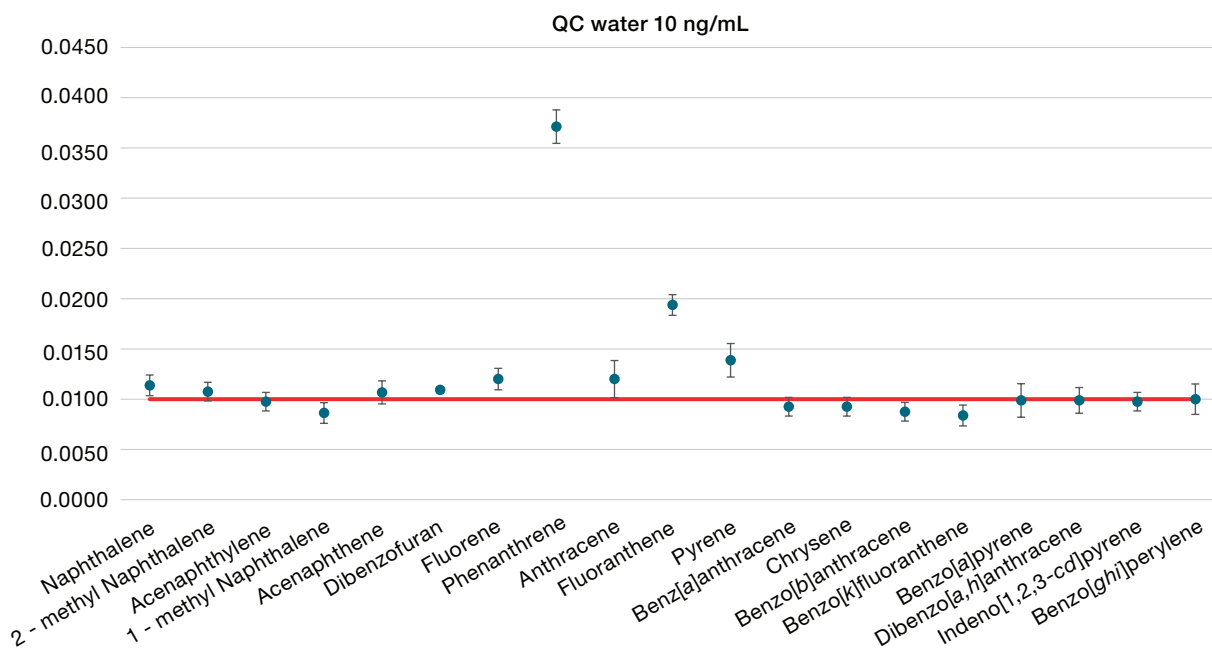


Figure A2. Average values of the calculated concentration of each compound in QC water at 10 ng/mL, with 2xStd error bars over the batch. Accuracy for some compounds, like phenanthrene, anthracene, fluoranthene and pyrene is not satisfactory, because QC was prepared from spiked matrix, which contained some analytes. Despite this bias, %RSD for all the compounds was <10%.

Table A3. Calculated concentrations of QC soil at 1,000 ng/mL during the robustness study, accuracy, standard deviation, and %RSD

QC soil 1,000 ng/mL	Calculated value	Target	Accuracy	2*STDEV	%RSD
Naphthalene	0.9405 0.9491 0.9452 0.9466 0.9588 0.9492 0.9476 0.939 0.9470	1.0000	95	0.01215	0.64
2 - methyl Naphthalene	0.9516 0.9691 0.9589 0.9522 0.9752 0.9705 0.9758 0.9487 0.9628	1.0000	96	0.02234	1.16
1 - methyl Naphthalene	0.9199 0.9443 0.937 0.9181 0.937 0.9247 0.9294 0.9068 0.9272	1.0000	93	0.02447	1.32
Acenaphthylene	0.902 0.8923 0.8726 0.8726 0.8973 0.9159 0.8576 0.8863 0.8871	1.0000	89	0.03760	2.12
Acenaphthene	0.9071 0.8848 0.9018 0.8967 0.8846 0.9094 0.893 0.9064 0.8980	1.0000	90	0.01969	1.10
Dibenzofuran	0.9277 0.9129 0.9249 0.9234 0.9133 0.9364 0.9097 0.9291 0.9222	1.0000	92	0.01867	1.01
Fluorene	0.9022 0.9161 0.8829 0.872 0.9315 0.8795 0.8375 0.9219 0.8930	1.0000	89	0.06207	3.48
Phenanthrene	0.9536 0.9431 0.9748 0.948 0.9534 0.9441 0.9585 0.9432 0.9523	1.0000	95	0.02139	1.12
Anthracene	0.9362 0.9222 0.9634 0.927 0.9382 0.93 0.9336 0.9111 0.9327	1.0000	93	0.03026	1.62
Fluoranthene	0.9268 0.9503 0.9599 0.9621 0.9497 0.9275 0.946 0.9535 0.9470	1.0000	95	0.02666	1.41
Pyrene	0.8226 0.8502 0.8495 0.8248 0.8869 0.9244 0.9216 0.8751 0.8694	1.0000	87	0.07940	4.57
Benz[a] anthracene	0.9093 0.9268 0.9513 0.927 0.932 0.9587 0.9517 0.9433 0.9375	1.0000	94	0.03317	1.77
Chrysene	0.9087 0.9298 0.927 0.9079 0.9158 0.921 0.9324 0.9169 0.9199	1.0000	92	0.01853	1.01
Benzo[b] fluoranthene	0.9014 0.9339 0.9128 0.9295 0.9596 0.9567 0.9849 0.9257 0.9381	1.0000	94	0.05467	2.91
Benzo[k] fluoranthene	0.9112 0.9501 0.8759 0.8538 0.9623 0.9754 0.9788 0.923 0.9288	1.0000	93	0.09263	4.99
Benzo[a]pyrene	0.9024 0.8974 0.9001 0.9019 0.9178 0.9072 0.9424 0.9144 0.9105	1.0000	91	0.02944	1.62
Dibenzo[a,h] anthracene	0.8925 0.8763 0.9141 0.891 0.8867 0.8564 0.8789 0.9064 0.8878	1.0000	89	0.03600	2.03
Indeno[1,2,3-cd] pyrene	0.8685 0.8495 0.8878 0.8773 0.8496 0.8228 0.8481 0.9008 0.8631	1.0000	86	0.05059	2.93
Benzo[ghi] perylene	0.8668 0.8443 0.8747 0.8595 0.8431 0.9412 0.8496 1.0073 0.8858	1.0000	89	0.11677	6.59

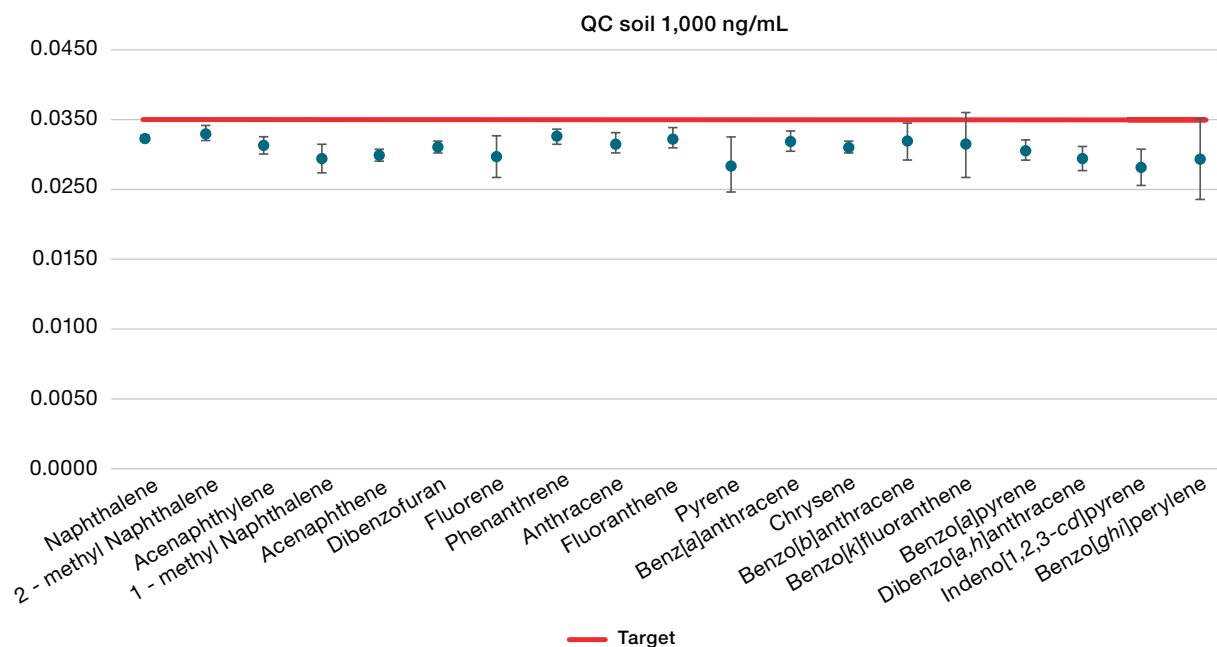


Figure A3. Average values of the calculated concentration of each compound in QC soil at 1,000 ng/mL, with 2xStd error bars over the batch

 Learn more at [thermofisher.com](https://www.thermofisher.com)

**For Research Use Only. Not for use in diagnostic procedures.** © 2021 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries. This information is presented as an example of the capabilities of Thermo Fisher Scientific products. It is not intended to encourage use of these products in any manners that might infringe the intellectual property rights of others. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representatives for details. **AN000455-EN 1221C**