

Application News

Gas Chromatograph HS-20 NX / Nexis™ GC-2060

Residual Solvent Analysis in Pharmaceuticals with N₂ Carrier Gas Using Nexis GC-2060 and HS-20 NX

— JP 18 and USP <467> / Water-Soluble Samples —

Shota Hayakawa, Toyohito Wada

User Benefits

- ◆ The newly designed FID delivers outstanding sensitivity, ensuring highly accurate and reliable analysis.
- ◆ Achieve full compliance with JP and USP guidelines using N₂ carrier gas, which is highly affordable and readily available.
- ◆ Easily analyzes *tert*-butyl alcohol and Cyclopentyl methyl ether, the Class 2 solvents added in the ICH Q3C (R8) guideline.

■ Introduction

For the analysis of residual solvents in pharmaceuticals, headspace gas chromatography (HS-GC) is widely used to target Class 1 and Class 2 solvents, in accordance with USP General Chapter <467> and JP18 (Japanese Pharmacopoeia, 18th Edition). These regulatory tests demand high sensitivity and reproducibility. The newly developed FID for the Nexis GC-2060 offers superior detection sensitivity compared to conventional models, enabling highly sensitive analysis. Additionally, due to recent helium supply shortages and rising costs, there is a growing demand for alternative analysis using nitrogen (N₂) as a carrier gas.

This application note presents the analysis of Class 1 and Class 2 water-soluble samples using the Nexis GC-2060 with N₂ carrier gas, compliant with Supplement II to the Japanese Pharmacopoeia. DMSO was used as the sample solvent.

* Some modifications were made to the USP and JP sample preparation protocols.

■ Analytical Conditions

The instrument configuration and analytical conditions for water-soluble samples are summarized in Table 1. To comply with USP and JP18 guidelines, Class 1 and Class 2 solvents were tested under two different methods (Procedure A and Procedure B) each utilizing distinct column types, column temperatures, and split ratios.

Table 1 Analytical Conditions for Water-soluble samples

GC Conditions (Procedure A / Procedure B)	
Model	: Nexis GC-2060
Detector	: FID (Flame Ionization Detector)
Column	: A) SH-I-624Sil MS (P/N 227-36066-02) (0.53 mm I.D. × 30 m, d.f.= 3.0 μm) B) SH-PolarWax (P/N 221-75972-30) (0.32 mm I.D. × 30 m, d.f.= 0.25 μm)
Column Temperature	: A) 40°C (20 min) – 10°C/min – 240°C (20 min) Total 60 min B) 50°C (20 min) – 6°C/min – 165°C (20 min) Total 59.17 min
Injection Mode	: A) Split 1:5 B) Split 1:10
Carrier Gas Controller	: Linear velocity (N ₂)
Linear Velocity	: 35 cm/sec
Detector Temperature	: 250°C
FID H ₂ Flow Rate	: 40 mL/min
FID Make up Flow Rate	: 30 mL/min (N ₂)
FID Air Flow Rate	: 170 mL/min

HS Conditions (Common to Procedure A and Procedure B)

Model	: HS-20 NX
Oven Temperature	: 80°C
Sample Line Temperature	: 110°C
Transfer Line Temperature	: 120°C
Vial Shaking Level	: Off
Vial Volume	: 20 mL
Vial Equilibrating Time	: 60 min
Vial Pressurizing Time	: 1 min
Vial Pressure	: 75 kPa
Loading Time	: 0.5 min
Load Equilib. time	: 0 min
Needle Flush Time	: 5 min
Injection Volume	: 1 mL

■ Analysis of Class 1 Standard Solution

Chromatograms obtained for the Class 1 standard solution using Procedure A and Procedure B are shown in Fig. 1 and 2, respectively. In addition, the signal-to-noise (S/N) ratio for each peak and the repeatability over six consecutive injections are summarized in Table 2 (for Procedure A) and Table 3 (for Procedure B).

In Procedure A, carbon tetrachloride showed an excellent S/N ratio of 23, with a relative standard deviation (%RSD) of 2% or less for individual peak areas. In Procedure B, 1,2-dichloroethane achieved an outstanding S/N ratio of 198, with a %RSD of 3% or less for individual peak areas, demonstrating highly favorable results.

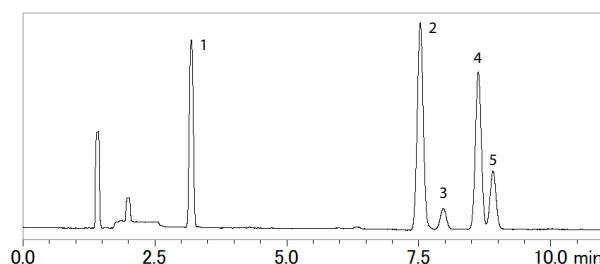


Fig. 1 Chromatogram of Class 1 Standard Solution (Procedure A)

Table 2 S/N Ratio and Repeatability of Class 1 Standard Solution (Procedure A)

Peak No.	Compounds	S/N ratio ave. ^{*1} (n = 6)	%RSD ^{*1} (n = 6)
1	1,1-Dichloroethane	331	1.23
2	1,1,1-Trichloroethane	339	1.26
3	Carbon tetrachloride	23	1.60
4	Benzene	195	1.94
5	1,2-Dichloroethane	67	0.91

*1 S/N ratios and %RSD are reference values and are not guaranteed.

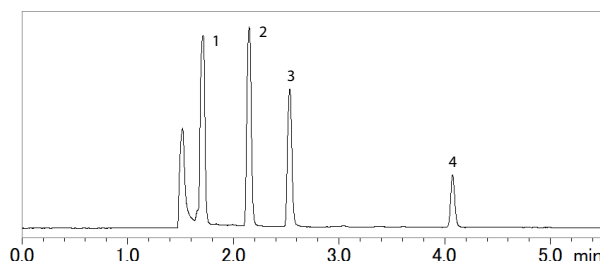


Fig. 2 Chromatogram of Class 1 Standard Solution (Procedure B)

Table 3 S/N Ratio and Repeatability of Class 1 Standard Solution (Procedure B)

Peak No.	Compounds	S/N ratio ave. ^{*2} (n = 6)	%RSD ^{*2} (n = 6)
1	1,1-Dichloroethane	157	1.73
2	1,1,1-Trichloroethane + Carbon tetrachloride	362	2.03
3	Benzene	389	1.87
4	1,2-Dichloroethane	198	1.15

*2 S/N ratios and %RSD are reference values and are not guaranteed.

■ Analysis of Class 2 Standard Solution

Chromatograms obtained using Procedure A and Procedure B are presented in Figure 3 and Figure 4, respectively (Class 2A: black, Class 2B: pink, TBA/CPME/MiBK: blue). Regarding system suitability, both methods yielded excellent results: Procedure A achieved a resolution of 1.0 or greater between the acetonitrile and dichloromethane peaks (Figure 3), while Procedure B achieved a resolution of 1.0 or greater between the acetonitrile and *cis*-1,2-dichloroethene peaks.



Note that the mixed standard sample containing *tert*-butyl alcohol (TBA), Cyclopentyl methyl ether (CPME), and Methyl isobutyl ketone (MiBK) was prepared separately to meet the specified concentrations.

■ Conclusion

Using the Nexis GC-2060 with nitrogen (N₂) carrier gas, we successfully achieved the analytical precision required by the Japanese Pharmacopoeia (JP18) and USP General Chapter <467>. The HS-GC method delivered excellent S/N ratios and high repeatability for the analysis of residual solvents in water-soluble pharmaceutical samples.

Related Applications

1. Residual Solvent Analysis in Pharmaceuticals with N₂ Carrier Gas Using Nexis GC-2060 and HS-20 NX — JP 18 and USP <467> / Water-Insoluble Samples — [Application News No. 01-01172-EN](#)

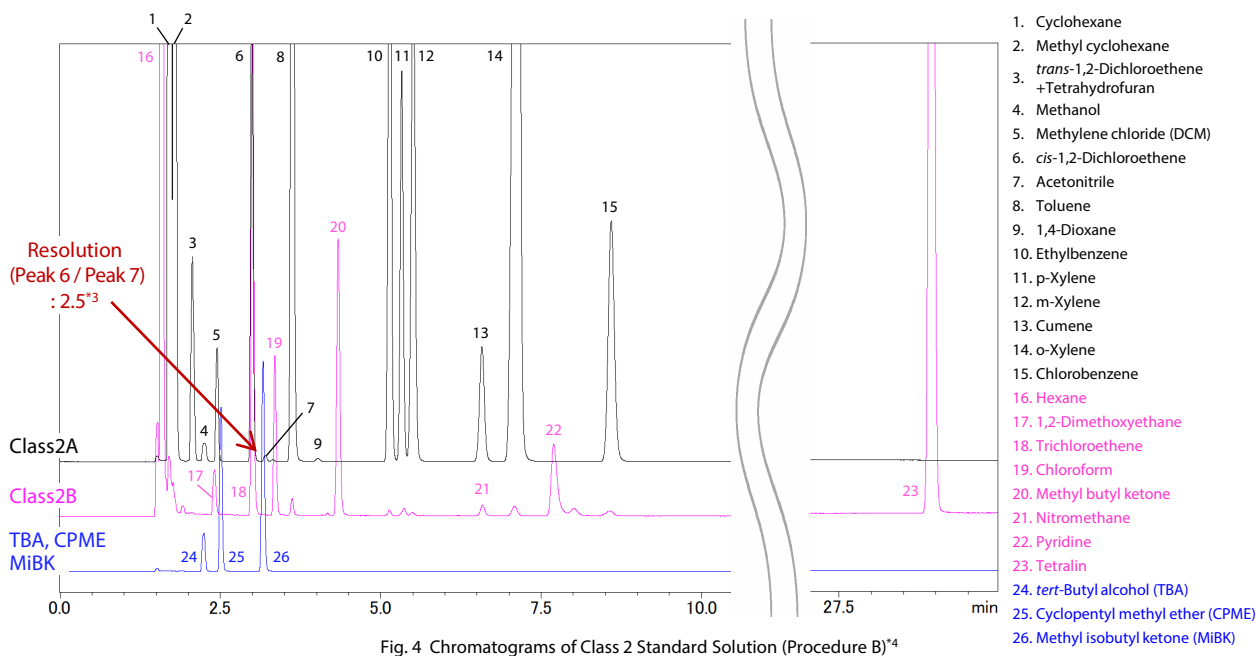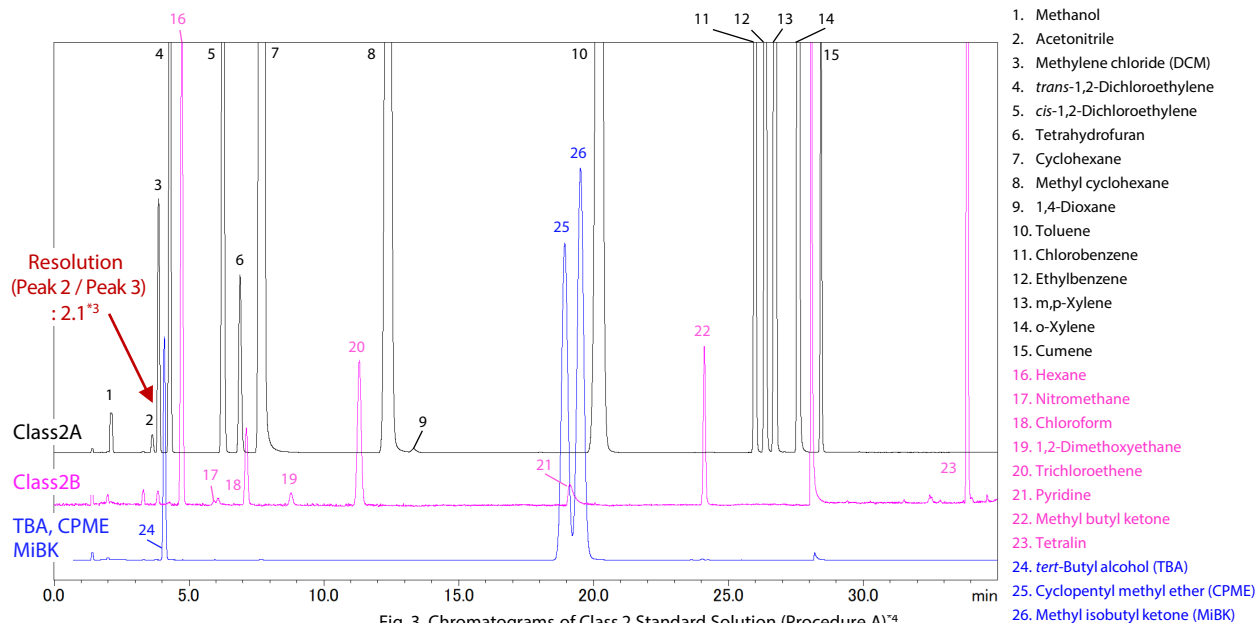

*³ The resolution values are reference values and are not guaranteed.

*⁴ The chromatograms in Figures 3 and 4 are displayed with different magnification scales.

Nexis is a trademark of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

01-01173-EN

First Edition: Aug. 2026



Shimadzu Corporation

www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. See <https://www.shimadzu.com/about/trademarks/index.html> for details.

Third party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

© Shimadzu Corporation, 2026

➤ Please fill out the survey

Related Products

Some products may be updated to newer models.



➤ Nexis GC-2060
Gas Chromatograph



➤ HS-20 NX series
Headspace Sampler

Related Solutions

➤ Pharmaceutical and
Biopharmaceutical

➤ Small Molecule
Pharmaceutical

➤ Life Science

➤ Price Inquiry

➤ Product Inquiry

➤ Technical Service /
Support Inquiry

➤ Other Inquiry