**CoreFocus**

P/N : 227-35032-01

Inlet Liner Removal Tool

**INSTRUCTION MANUAL**

The Inlet Liner Removal Tool is designed to simplify and efficiently facilitate the removal of gas chromatography (GC) inlet liners.

Applicable liner outer diameter : 5 mm

Recommended operating temperature : Room temperature


**! CAUTION**

Please use this tool only after SPL has cooled down to room temperature.

There is a risk of burns when SPL is at high temperatures. Additionally, using this tool at high temperatures may accelerate its deterioration.

**【The procedure for removing the liner】**

1. While holding down the injection port, turn the lever counterclockwise.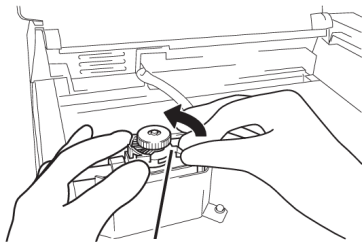

Lever


2. Slowly lift the injection port.

**Note**

Please lift the injection port slowly and directly upwards with your hand to prevent damage to the glass liner.

**Tips**

If the O-ring is stuck on the injection port, please use a pair of tweezers to carefully remove the O-ring before lifting the injection port to prevent any damage to the glass liner.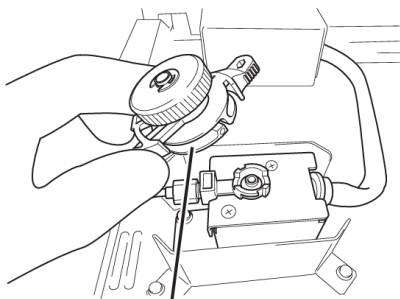


Injection Port

3. Insert the tool into the liner that is protruding from the SPL.


**Note**

Please be careful not to damage the sealing surface of the O-ring on the sample injection port.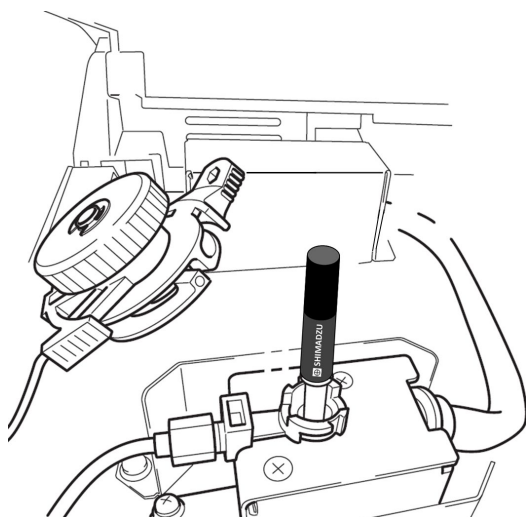

**Tips**


If the O-ring is stuck due to heat, please use a pair of tweezers to carefully tear the O-ring away from the main body via the two grooves on both sides.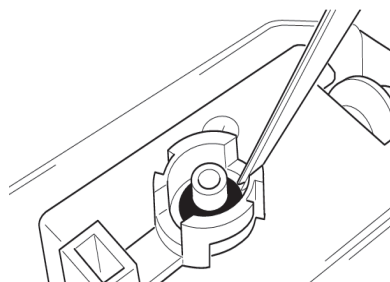

4. Pull it straight up to remove it.

**Note**

There is a possibility of the liner breaking if you pull it out diagonally.

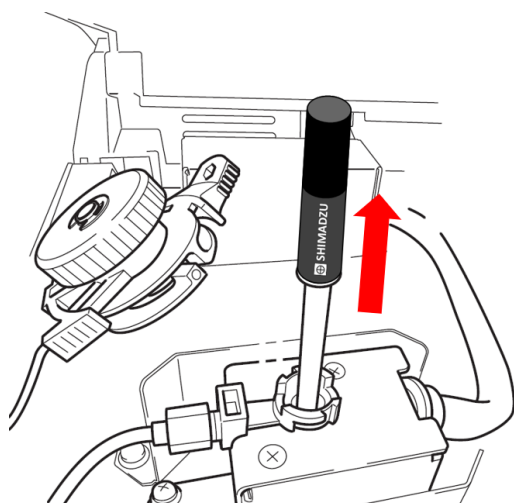5. Remove the tool from the extracted liner.

**Note**

Please be careful as the liner may be hot.

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