

Biotage® Horizon 5000

Release Notes



What's New in Biotage® Horizon 5000 1.1.0.6

Enhancements

- » **Solvent Dispensing Calibration.** The “Density” field on the Solvent Library Screen is renamed to “Density Factor” to reflect the users’ ability to use values other than the standard reagent density. Historically, the field is used to control the amount of reagent delivered for conditioning and elution in a method. Nominally set to the density (in g/mL) of the reagent being used, this value can now be customized to allow for more precise amounts of reagent to be dispensed depending on factors such as the reagent or reagent mixture and the ambient laboratory conditions.
- » **Additional Error Reporting.** If the PDF reader that is set up is not found by the software on the PC, a message box is now displayed stating: “.pdf reader not found”.

Bug Fixes

- » **Corrected Communication Errors** which typically result in situations where a station that is asked to start, does not. Previously, to compensate for this error, the user would need to add the sample to the next run or restart all stations being used, ensuring that all requested stations have started correctly.

Compatibility Information

- » It is important with the recent 5000 software release v1.1.0.6, to update the USB driver and the firmware in the 5000 extractors to v1.0.0.3 to prevent communication issues.

For more information, please contact your local Biotage representative. See contact information on our website www.biotage.com.

EUROPE

Main Office: +46 18 565900
Toll Free: +800 18 565710
Fax: +46 18 591922
Order Tel: +46 18 565710
Order Fax: +46 18 565705
order@biotage.com
Support Tel: +46 18 56 59 11
Support Fax: + 46 18 56 57 11
eu-1-pointsupport@biotage.com

NORTH & LATIN AMERICA

Main Office: +1 704 654 4900
Toll Free: +1 800 446 4752
Fax: +1 704 654 4917
Order Tel: +1 704 654 4900
Order Fax: +1 434 296 8217
ordermailbox@biotage.com
Support Tel: +1 800 446 4752
Outside US: +1 704 654 4900
us-1-pointsupport@biotage.com

JAPAN

Tel: +81 3 5627 3123
Fax: +81 3 5627 3121
jp_order@biotage.com
jp-1-pointsupport@biotage.com

CHINA

Tel: +86 21 68162810
Fax: +86 21 68162829
cn_order@biotage.com
cn-1-pointsupport@biotage.com

KOREA

Tel: +82 31 706 8500
Fax: +82 31 706 8510
korea_info@biotage.com
kr-1-pointsupport@biotage.com

INDIA

Tel: +91 22 4005 3712
india@biotage.com

Distributors in other regions are listed on www.biotage.com

Literature Number: PPS594

© 2019 Biotage. All rights reserved. No material may be reproduced or published without the written permission of Biotage. Information in this document is subject to change without notice and does not represent any commitment from Biotage. E&OE. A list of all trademarks owned by Biotage AB is available at www.biotage.com/legal. Other product and company names mentioned herein may be trademarks or registered trademarks and/or service marks of their respective owners, and are used only for explanation and to the owners’ benefit, without intent to infringe.